

Designing a Dry Material Feeder

CHAD M. LORENSEN, Pharmaceutical Industry Business Development Manager, & ANDREAS LUDWIG, Mechanical Engineering Manager, Schenck AccuRate Inc.



Metering dry powders in the pharmaceutical industry is not a new idea. Several common applications have been utilizing gravimetric and volumetric dry material feeders for many years. Until recently, these feeders were customized versions of industrial or food grade equipment or one-off type machines engineered for a specific application. Driven by increasing demand, an industry leading equipment manufacturer embarked on an effort to create a standard dry material feeder engineered to meet the pharmaceutical industries' needs.

As part of this process, common pharmaceutical applications utilizing feeders were identified: milling/particle size reduction; mix tank and reactor charging; mixing, pharmaceutical grade plastics extrusion; tablet coating; inter-process packaging; and continuous granulation. Once the applications were identified, a multi-functional research and design team interviewed technical people from companies who built equipment specifically for the pharmaceutical industry as well as companies who would use the equipment in their own production efforts. This research was conducted in North America as well as Europe to provide a global picture of the requirements for the product.

Many design hurdles were discovered during the interviews and review of feeding equipment requirements for these applications. The hurdles were made even more complicated by the fact that there were no widely accepted industry standards for the manufacture of processing equipment. Each manufacturer and end user had

Designing a Dry Material Feeder

Published on Chem.Info (<http://www.chem.info>)

different standards they felt were important.

The major criteria addressed during design included: equipment capacities and rates; FDA approved materials of construction, surface finishes of welds and metals used in fabrication; cleaning processes used on the equipment during production including frequent assembly and disassembly; validation; and feeding accuracy. The following paragraphs elaborate on the various design criteria for the equipment.

Equipment Capacities & Rates

During interviews, the research and design team studied capacities of typical downstream processing equipment that would be paired with the dry material feeder. Based on the most common sizes of these machines, two distinct application feed rate ranges were determined: 20 - 2,000 g/hr and 0.5 - 150 kg/hr. These ranges were used to determine the appropriate size and design of equipment. For the purpose of this paper, we will review the design process for the machine at the higher range of 0.5 - 150 kg/hr.



FDA-Approved Materials of Construction

The interviews determined 316 stainless steel is a preferred material for equipment in the pharmaceutical industry because of its corrosion, chemical and thermal resistance. EPDM was found to be a widely accepted polymer used for components that have to be elastic, non porous and FDA compliant. The design team worked with various suppliers and molders of EPDM to adapt the polymer for use in the flexible hopper. Flexibility and temperature resistance made it ideal for this application. Additionally, the cost effective nature of the flexible hopper allows it to be used as a disposable component versus the higher cost of a typical cleaning operation.

Designing a Dry Material Feeder

Published on Chem.Info (<http://www.chem.info>)

Other non-metal materials of construction include PTFE, PET-P and Nylon.

Even though the proper selection of construction materials is essential, geometry and surface finish are no less important when designing a pharmaceutical grade machine.

Welding, Surface Finishes & General Design

Existing experience from food and dairy feeding applications dictated that all welds must be continuous, non porous, and without cracks or undercuts to ensure the machine can be easily cleaned. Utilizing larger surfaces versus complex designs with undercuts give the machine easy access for cleaning and



maintenance. Additionally, larger, smooth surfaces help eliminate the potential for water or solvent to remain in crevices.

Studies indicated 32 microinch roughness average (Ra) was a widely accepted surface finish however; certain applications dictated finishes up to 16 microinch Ra. Consequently, standard configurations for 32 Ra and 16 Ra were developed.

If complex design cannot be avoided, then providing easy disassembly for cleaning purposes must be strongly considered. Examples include the use of stainless steel tri-clover type clamps to secure the extension hopper and discharge nozzle as well as a continuously welded housing, agitation arm and other peripheral components.

Cleaning Processes Used on the Equipment

Since product contamination from foreign particles must be strictly avoided, materials of construction have to withstand aggressive cleaning solutions and high temperatures to handle regular maintenance without deterioration or corrosion while maintaining FDA compliance.

Common sense dictates fewer components in the product contact zone will reduce cleaning efforts. External agitation of the material is a good example of this concept and was selected by the design team over internal agitation.

Designing a Dry Material Feeder

Published on Chem.Info (<http://www.chem.info>)



Validation

Since pharmaceutical companies are required to verify the product contact materials, it is essential for the manufacturer to keep accurate records of raw material certifications. The manufacturer implemented procedures to procure raw materials with certification and track them through their plant and manufacturing process to guarantee proper material composition. A standard process was created to generate IQ/OQ documentation for each individual feeder manufactured.

Feeding Accuracy

Since pharmaceutical grade feeders are often used to feed active pharmaceutical ingredients into a continuous process, feeding accuracy is essential.

Many of those ingredients have different flow behaviors including cohesive and adhesive properties making variable feeder settings advantageous.

The design team created an agitation system enabling variable speed and penetration depth. These features allow the user to find the ideal setup for an optimal product flow into the discharge element. Consistent helix filling is critical for good discharge behavior, resulting in precise accuracies. The control algorithm in gravimetric feeders allows for continual helix speed adjustments depending on the loss in weight over time which again increases the feeding accuracy. The manufacturer's proven gravimetric control system was utilized to facilitate this functionality.

Designing a Dry Material Feeder

Published on Chem.Info (<http://www.chem.info>)



When feeding cohesive materials like those typically found in the pharmaceutical industry, often result in an uneven discharge behavior. The material being fed drops off in lumps instead of creating an even flow. One way to prevent this can be a screened discharge nozzle, which breaks up agglomerations of bulk materials and produces a more uniform product flow.

The design team gave special considerations for the design of a screened discharge nozzle that was used to facilitate disassembly and cleaning.

When designing or specifying a pharmaceutical feeder or feeding system, it is important to consider basic design principles, materials being fed, cleaning processes used, and FDA requirements. Implementing a properly designed feeder into your process can save time and reduce overall manufacturing costs.

For more information, please e-mail mktg@accuratefeeders.com [1] or visit www.accuratefeeders.com [2].

Source URL (retrieved on 04/19/2015 - 10:45am):

<http://www.chem.info/articles/2010/09/designing-dry-material-feeder>

Links:

[1] <mailto:mktg@accuratefeeders.com>

[2] <http://www.accuratefeeders.com/>